



December 2021

Xsens

Release notes MVN
2021.2

MVN 2021.2 release notes

New versions	MVN: 2021.2 (Dec 2021) Awinda firmware: (unchanged) AW-A2 v4.6.0, AW-DNG2(-ANT) v4.6.0, MTw2 v4.6.0 Link firmware: (unchanged) BPACKv1 v1.6.1, BPACKv2 v1.6.1, MTx2 v1.6.1 FWU: 4.3.1
Previous versions	MVN: 2021.0.1 (Jun 2021) Awinda firmware: AW-A2 v4.4.0, AW-DNG2 v4.4.0, MTw2 v4.4.0 Link firmware: BPACKv1 v1.6.0, BPACKv2 v1.6.0 MTx2 v1.6.0 FWU: 4.3.1
Release date	December 2021
Operating system	Windows 10, 64-bit
Meaning of icons	+added, ^changed, !note, #fixed, -removed

1 Introduction

This document describes the new features, improvements, and bug fixes in the new release 2021.2 for Xsens software: MVN Record, MVN Analyze Plus, MVN Analyze Pro, MVN Animate Plus and MVN Animate Pro.

2 Major improvements and new features

2.1 MVN

- + Easier and faster On-the-Go calibration
- + Up to 50% faster HD reprocessing
- + EMG data added to C3D exporter
- + Improved workflow for OBR import
- + Improved MotionCloud integration and workflow
- + MVN Basic renamed to MVN Record
- + Xsens gloves by Manus enabled for MVN Record
- + MVN Animate- and Analyze Plus now allows to connect 2 systems simultaneously

3 Bug fixes

3.1 MVN

- # Resolved command-line hanging when trying to reprocess a read-only file
- # Resolved command-line export to MVNX inserting frames
- # Resolved MVN crashing when swapping trackers
- # Resolved MVN crashing when playing during reprocessing
- # Avatar's name in viewport is now displayed correctly in Japanese & Korean
- # MOBU can receive object tracking data when MVN is using non-English language
- # Calibration is saved when a suit gets disconnected
- # Limited file name preventing MVN from crashing

- # FBX exporter output preset fix
- # Changes in a mvns file can be saved correctly
- # Viewport fix when pressing Alt+ set to Free Move
- # Resolved "Stop Network Streaming to target"

4 Minor updates

4.1 MVN

- + Improved Network Streamer
- + Ability to use Vive Z-value in no level and multilevel (non live only)
- + Improved R(OBR) body dimension configuration window
- + Install plugins is now optional in the MVN installer
- + Accept different formats for time code time sync

5 Known Issues

None